

## Acid Phenol:Chloroform MB Grade

Store at 4°C or -20°C.

<b>Catalog #:</b>	AM9720
<b>Amount:</b>	100 mL
<b>Product Description:</b>	Acid phenol:chloroform, premixed with isoamyl alcohol (125:24:1 phenol:chloroform:isoamyl alcohol).
<b>Appearance:</b>	Clear colorless liquid
<b>Chemical Analysis:</b>	Purity >99.0% pH 4.3–4.7
<b>Caution:</b>	<b>Phenol is very corrosive and will severely burn the skin. Safety precautions such as gloves, protective eyewear, a lab coat, and working in a fume hood are critical.</b>
<b>Storage Conditions:</b>	4°C or -20°C. Store in amber or foil-wrapped bottles, and do not open frequently, to avoid oxidation and breakdown products. To prolong the shelf life, store aliquots in 50 mL tubes at -20°C.

---

### USER INFORMATION

---

#### General Information:

Phenol is a crystalline solid at room temperature. It is normally either saturated with an aqueous buffer or dissolved in chloroform to make it liquid. Phenol extraction is a commonly used method for deproteinization of nucleic acids [1, 2, 3]. Most proteins are more soluble in phenol than in the aqueous phase. Conversely, nucleic acids are more soluble in the aqueous phase. Centrifugation of the mixture will yield two phases; the lower phase is the organic phase and will contain the protein, usually as a white flocculent at the interface. The upper aqueous phase will contain nucleic acids.

Phase partitioning of nucleic acids is pH dependent [4, 5]. At pH 4–6 DNA will be retained in the organic phase and interface, leaving the RNA in the aqueous phase. Isolation of RNA [5, 6] from biological material is often done with an acid-phenol [6, 7]. For DNA isolation, a pH of 7.5–8.0 is required, and both DNA and RNA will partition into the upper aqueous phase. The pH dependence of DNA phase partitioning makes it necessary to raise the pH of many in vitro procedures prior to phenol extraction. For applications which require an alkaline pH, such as DNA isolation, use phenol buffered to pH ~7.9.

Salt concentration also has an effect on the efficiency of phenol extraction. RNA has traditionally been extracted at high salt concentration (0.2–0.5 M NaCl) to reduce the effect of endogenous ribonucleases. With the use of more effective ribonuclease inhibitors, such as guanidinium salts, optimal RNA yields are obtained in an extraction buffer with low or no salt present.

Chloroform is mixed with phenol to increase the efficiency of nucleic acid extractions. The increased efficiency of extraction is due to chloroform's ability to denature proteins, thus aiding separation of nucleic acid from protein. Phase separation of the extracted solution is also enhanced, thereby assisting in the removal of the aqueous phase with minimal cross contamination from the organic phase. Addition of chloroform to the phenol also aids in the removal of lipids. Often isoamyl alcohol (IAA) is added to the phenol:chloroform to prevent foaming.

#### Applications:

##### Notes on Use of Phenol

**To measure the pH of phenol:** Measurement of phenols with pH paper is not accurate due to breakdown of indicator chemicals on the paper. Mix 2 mL of the organic phase with 8 mL of methanol and 10 mL of water. Measure the pH of the entire sample with a pH meter.

Preparation of crystalline phenol for use is a time-consuming and often hazardous procedure. The Ambion® series of Saturated Phenols eliminates these handling problems.

Product	Catalog #	Applications
Saturated Phenol pH 6.6/7.9 ± 0.2	AM9710 – 100 mL AM9712 – 400 mL	Extraction of nucleic acids Add alkaline buffer (included) for pH 7.9
Acid Phenol:Chloroform pH 4.7	AM9720 – 100 mL AM9722 – 400 mL	RNA isolation Removing DNA from in vitro transcription reactions
Phenol:Chloroform:IAA pH 6.6/7.9 ± 0.2	AM9730 – 100 mL AM9732 – 400 mL	Extraction of nucleic acid, when phenol:chloroform is desired. Add alkaline buffer (included) for pH 7.9

**References:**

1. Ausubel FM, Brent R, Kingston RE, Moore DD, Seidman JG, Smith JA, Struhl K, editors. (1987) *Current Protocols in Molecular Biology*, p 2.1.3–2.1.4, 15.3.8–15.3.11.
2. Wallace DM. (1987) Large- and small-scale phenol extractions. *Methods in Enzymology* **152**: 33–41.
3. Sambrook J, Fritsch EF, and Maniatis T. (1989) *Molecular Cloning: A Laboratory Manual*, 2nd Ed.
4. Perry RP, La Torre J, Kelley DE, Greenberg JR. (1972) On the lability of poly(A) sequences during extraction of messenger RNA from polyribosomes. *Biochim. Biophys. Acta* **262**: 220–226.
5. Brawerman G, Mendecki J, Lee SY. (1972) A procedure for the isolation of mammalian messenger ribonucleic acid. *Biochemistry* **11**: 637–641.
6. Chirgwin JM, Przybyla AE, MacDonald RJ, Rutter WJ. (1979) Isolation of biologically active ribonucleic acid from sources enriched in ribonuclease. *Biochemistry* **18**:5294–5299.
7. Chomczynski P and Sacchi N. (1987) Single-step method of RNA isolation by acid guanidinium thiocyanate-phenol-chloroform extraction. *Anal Biochem* **162**:156–159.

---

**QUALITY CONTROL**

---

**Functional Testing:**

Tested for phase separation when combined with an aqueous solution.

The pH of Acid Phenol:Chloroform in a 10% methanol solution is 4.4 ± 0.2.

---

**OTHER INFORMATION**

---

**Material Safety Data Sheets:**

Material Safety Data Sheets (MSDSs) can be printed or downloaded from product-specific links on our website at the following address: [www.ambion.com/techlib/msds](http://www.ambion.com/techlib/msds). Alternatively, e-mail your request to [MSDS\\_Inquiry\\_CCRM@appliedbiosystems.com](mailto:MSDS_Inquiry_CCRM@appliedbiosystems.com). Specify the catalog or part number(s) of the product(s), and we will e-mail the associated MSDSs unless you specify a preference for fax delivery. For customers without access to the internet or fax, our technical service department can fulfill MSDS requests placed by telephone or postal mail. (Requests for postal delivery require 1–2 weeks for processing.)

**Warranty and Liability:**

**For research use only. Not for use in diagnostic procedures.**

Applied Biosystems is committed to delivering superior product quality and performance, supported by industry-leading global service and technical support teams. Warranty information for the accompanying consumable product is available at [www.ambion.com/info/warranty](http://www.ambion.com/info/warranty) in "Limited Warranty for Consumables," which is subject to the exclusions, conditions, exceptions, and limitations set forth under the caption "EXCLUSIONS, CONDITIONS, EXCEPTIONS, AND LIMITATIONS" in the full warranty statement. Please contact Applied Biosystems if you have any questions about our warranties or would like information about post-warranty support.

Information in this document is subject to change without notice. Applied Biosystems assumes no responsibility for any errors that may appear in this document.

Applied Biosystems disclaims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall Applied Biosystems be liable, whether in contract, tort, warranty, or under any statute or on any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

**Trademarks, Patents, and Licensing:**

Applied Biosystems, AB (Design), and Ambion are registered trademarks of Applera Corporation or its subsidiaries in the US and/or certain other countries. All other trademarks are the sole property of their respective owners.

© 2008 Ambion, Inc. All rights reserved. 4386643A